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Appearance	Colorless to slightly yellow Liquid/Suspension
Physical Form	Colloidal solution of silica coated with polyvinylpyrrolidone (PVP)
Osmolality	< 25 mOsm/kg
Viscosity Max.	15 cP
Conductivity	<100 mS/m

pH Range	8.5–9.5
Density Max.	1.135 g/ml
Application	Used in balanced salt solutions, physiological saline or 0.25 M sucrose. Used for basic research applications. Used for density gradient centrifugation of cells, viruses and subcellular particles. Used for formation of gradients either by the use of convenient gradient mixers or by high speed centrifugation.
Estimated Shelf Life from Manufacture Date	5 years
Storage Conditions	4–30°C
Stability	Percoll is provided as solution which is stable for at least 5 yr when stored between 4°C and 30°C

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此公式中, V_x = volume of diluting medium (ml); V_0 = volume of undiluted Percoll (ml)

p_0 = density of Percoll (1.130+0.005g/ml);

p_{10} = density of 1.5M NaCl=1.058g/ml(minor differences for other salts);

density of 2.5M sucrose=1.316g/ml(minor differences for other additives)

ρ_i = density of SIP solution produced(g/ml)

Thus, for SIP in saline, $\rho_i=1.123\text{g/ml}$ and for SIP in sucrose, $\rho_i=1.149\text{g/ml}$, assuming $\rho_0=1.130\text{g/ml}$.

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$$V_y = V_i \frac{(\rho_i - \rho)}{(\rho - \rho_y)}$$

此公式中， V_y = volume of diluting medium in ml; V_i = volume of SIP in ml;

ρ_i = density of SIP in g/ml; ρ = density of diluted solution produced in g/ml;

ρ_y = density of diluting medium in g/ml

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